



PurMop BLACK®

F1-s

Cleanroom Mop –
Performance Line

Sterile | Fillpack (ready to
fill with fluid)

Polyester-Microfiber

Item no. 2112046

STERILE

ISO	3	4	5	6	7	8
GMP			A/B		C	D

The PurMop BLACK® F1-S is a fillpack with cleanroom mops of the latest generation. Inside are two mops of the D1 series. These convince with proven polyester microfiber cleaning performance, the Wings technology for better corner cleaning as well as a new size, which results in a higher area performance. The fillpack is a chemical-resistant pouch with a lid into which the fluid of choice is filled. A wide range of cleaning and disinfectant agents can be added, for which today there is no presaturated mop available. The pouch is resistant to common bactericides, sporicides and detergents used in cleanrooms. Together with the PurMop BLACK® carbon fiber mop systems, the F1-S provides an unprecedented level of ergonomics, performance and safety.

- Pouch with screw cap: ready for saturation with the fluid of your choice
- Pouch and mops are resistant to common cleanroom bactericides, sporicides and detergents
- “BLACK48” – new size suitable for PurMop BLACK® carbon fiber mop systems

Application ceiling / wall / floor

Industry sectors pharma / medical /
pharmacy / laboratory / mechanics

APPLICATION, USAGE

For single cleanroom application. Recommended for cleaning and disinfection of sterile GMP cleanrooms.

SIZE, WEIGHT, COLOR

48,0 x 12,0 cm (±0,5 cm) | 45 g (± 10 %) white, orange binding

FRONT MATERIAL

80% polyester microfiber /
20% polyester

BACKING MATERIAL

100% polyester nonwoven

EDGES

100% polyester overlock (bound)

FILLING QUANTITY

600 ml (for floors)
350 ml (for ceilings and walls)

SURFACE COVERAGE

Up to 40 m² (per fillpack–floors)
Up to 30 m² (per fillpack–ceilings/walls)

CHEMICAL COMPATIBILITY

Against acids: resistant
Against alkalis: resistant

ABRASION RESISTANCE

≥ 20.000 movements (9 kPa)

STERILITY

Gamma irradiated, SAL 10⁻⁶,
validated according to ISO 11137

SHELF LIFE

3 years

RE-USE, RECYCLING

Recyclable

STORAGE

General storage conditions apply, protect from harmful influences (fluids, sun rays or UV light, contamination, pressure), recommended storage temperature 5–40 °C

PACKING

2 mops / pack
45 packs / carton [90 mops]

CARTON SIZE (L/W/H)

CARTON WEIGHT
60 x 50 x 40 cm | 8,7 kg

PALLET

20 cartons / pallet
(120 x 100 x 223 cm)

CUSTOMS TARIFF NUMBER

COUNTRY OF ORIGIN
63071090,
Socialist Republic of Vietnam

COMPATIBLE ACCESSORIES

- PurMop BLACK® C1 & C2
- Carbon Fiber Mop Systems
- Funnel (Material: Polypropylene, Diameter: 120 mm, Item no. 2112060)
- Beaker (Volume: 1 l, Material: Polypropylene, Printed black scale, Item no. 2112059)



BLACK
PERFOR
MANCE
LINE



Hydroflex Group GmbH . Kiesacker 19 . 35418 Buseck . Germany

T +49 6408 50435-0 . F +49 6408 50435-20 . info@hydroflex-group.com . hydroflex-group.com

PurMop BLACK® F1-s

INSTRUCTIONS FOR USE

DESCRIPTION

Open the screw cap of the pouch and remove the sealing. Fill in the fluid and close the pouch again with the screw cap.

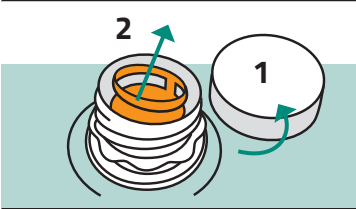
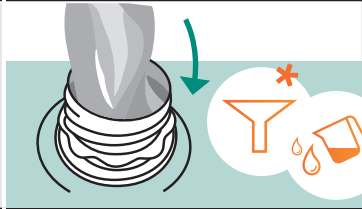
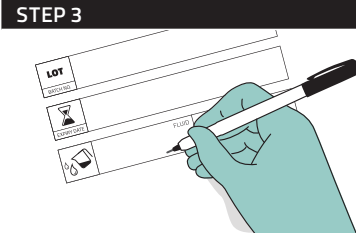
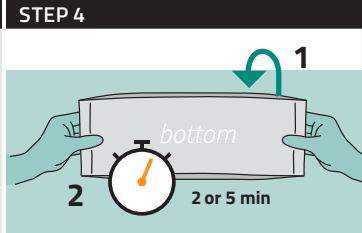
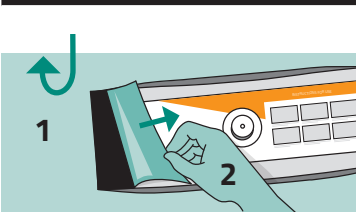
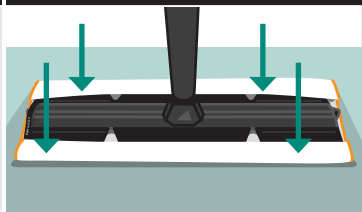
Fluid volume: 600 ml (for the floor) or 350 ml (for the ceiling and walls)

Note the filling date and the fluid used on the label.

For a complete saturation of the mops, turn the closed pouch upside down and wait 2 minutes for floors or 5 minutes for ceilings/walls.

After the waiting time, turn the pouch back and open it sideways on the left edge of the packaging by tearing open the peelable flap. The mops can be picked up directly from the fully opened packaging using the mop holder. The pouch cannot be resealed after opening the sealing seam.

Recommended: The compatibility of the chemical formulation of the fluid and the surface to be cleaned or disinfected should be checked before adding the liquid to the dry mops.

STEP 1	STEP 2
 - Unscrew the lid (1) - Remove sealing (2).	 - Add 600 ml of fluid (for floor) or 350 ml (for ceilings and walls) and close the lid.
STEP 3	STEP 4
 - Note the fluid used and the filling date on the label.	 - Turn the closed pouch upside down (1) and wait (2): 2 minutes for floors or 5 minutes for ceilings/walls.
STEP 5	STEP 6
 - Turn the pouch back to top (1) and open it by tearing from edge (2).	 - Take up mop with PurMop BLACK® Carbon Fiber Mop System.

* Funnel (Item No. 2112060) and Beaker (1 l, Item No. 2112059) available as optional accessories